

**How to get free academia license**

**for**

**TopSpin 3.6.X & CMC-assist 2.X**

2019.07.16

# Step 1: Register for Bruker users account.

<https://www.bruker.com/about-us/register.html>

Register - About Us | Bruker

Products Applications Service News Events About Us Login Contact

Home - About Us - Register

## Register for Bruker users account

Welcome to Bruker quick registration portal. To download Bruker software, you are required to have an account. If you don't have an account with us yet, you can start now by filling out the form below.

Please also note that the personal information that you enter in the form below may be screened against relevant sanction lists to ensure compliance with applicable export control laws.

\*= These are mandatory fields

First Name: \*

Last Name: \*

Country or Area: \* Taiwan

Email Address: \*

Create Password: \*

Confirm Password: \*

Password must be at least 6 characters long

I am interested in the following newsletters:

- Magnetic Resonance Academia and Research eNews
- Mass Spectrometry
- Atomic Force Microscopy Materials Journal Club
- Fluorescence Microscopy Journal Club
- Preclinical Imaging eNews
- Microbiology & Diagnostics
- Magnetic Resonance Education eNews
- Magnetic Resonance Pharma eNews
- Magnetic Resonance Clinical eNews
- Magnetic Resonance Food and Beverages eNews

By submitting this form I confirm that I would like to be contacted by Bruker. I have taken note of the Bruker Website [Privacy Notice](#) and I agree to the [Terms of Use](#) of the Bruker website. I understand that I may unsubscribe at any time.\*

I'm not a robot

reCAPTCHA  
Privacy - Terms

f  
t  
in  
e

## Step 2: Login using your account , and click “Free Academia license”.

<https://www.bruker.com/service/support-upgrades/software-downloads/nmr.html>

Home . Service . Support & Upgrades . Software Downloads . NMR

### NMR Software & Downloads

This section is available for customer only, registered users can access to NMR software download, upgrades and patches for Windows, Linux and macOS.

If you don't have an account, please [register for an account](#). If you already have an account please [log in here](#) to access the download section containing all listed software products below.

#### Software Downloads

If you are logged in, you can download the software below by clicking on the link:

- TopSpin for Windows, Linux and macOS  
*TopSpin installation includes following products:*
  - IconNMR
  - SmartDriveNMR
  - CMC-q
  - CMC-se
  - CMC assist
  - PotencyMR
  - FBS
  - TopSolids
  - BioTop

#### Generate License Code

TopSpin processing software is free for academia and government institutions. And you can get an Evaluation license for the full Bruker Suite [here](#) (requires login/ quick registration).

- [Free Academia license](#)
- [Free Evaluation license](#)

#### License Request

Request NMR processing and acquisition licenses

[Request Form](#)

Latest versions of Bruker NMR software use the CodeMeter licensing system. We offer a possibility of a complimentary exchange of existing FlexLM licenses for corresponding CodeMeter license tickets.

[NMR License Exchange Form](#)

1. Log in

2. Click “Free Academia license” .

## Step 3: Now, click "Request personal license ticket now" .

The screenshot shows a web browser window with the URL [https://www.bruker.com/service/support-upgrades/software-downloads/nmr/free-topspin-processing-...olicensesgeneration\[controller\]=DemoLicensesGeneration&cHash=3f31f0da1fc6e9f7244fc60490596194](https://www.bruker.com/service/support-upgrades/software-downloads/nmr/free-topspin-processing-...olicensesgeneration[controller]=DemoLicensesGeneration&cHash=3f31f0da1fc6e9f7244fc60490596194). The page features the Bruker logo and a navigation menu with links for Products, Applications, Service, News, Events, About Us, Logout, and Contact. A search bar is also present. The main content area is titled "Free TopSpin Processing for Academia" and includes a "Contact" button. Below this, a breadcrumb trail reads "Home . Service . Support & Upgrades . Software Downloads . NMR . NMR TopSpin License for Academia". The main heading is "NMR TopSpin License Processing for Academia", followed by a paragraph explaining the free access to TopSpin software for academic users. A prominent blue button labeled "Request personal license ticket now" is highlighted with a red arrow and the text "Click here". Below this button are three links: "Download for Windows", "Download for Linux", and "Download for Mac OS". A "Disclaimer" section is located at the bottom left, and social media icons for Facebook, Twitter, LinkedIn, and Email are on the bottom right.

Free TopSpin Processing for Academia

Contact

Home . Service . Support & Upgrades . Software Downloads . NMR . NMR TopSpin License for Academia

### NMR TopSpin License Processing for Academia

In order to provide students, researchers and teachers with unlimited access to the best tools for off-line NMR processing, Bruker is making their market leading NMR processing software TopSpin available free of charge for all academic users. In addition to Bruker NMR data, the software is able to process data acquired on Agilent, Varian and Jeol systems, as well as read any data provided in JCAMP format. We are happy to offer you a free Academia license for **TopSpin 4 and TopSpin 3.6** ("Academia license").

**Request personal license ticket now**

[Request personal license ticket now](#)

[Download for Windows](#)   [Download for Linux](#)   [Download for Mac OS](#)

#### Disclaimer

Access to Academia licenses is restricted to academic customers. By using this web page for license requests you confirm that you are employed by or studying at an academic- or governmental institution and that you use the software exclusively in this context. Your request and the license generation will be logged. If you do not agree with this, please do not continue.

In the next step a personal ticket code will be generated for you. With this ticket code you can activate the Academia license on a compatible computer of your choice and use the software:

[https://www.bruker.com/service/support-upgrades/software-downloads/nmr/free-topspin-processing-...olicensesgeneration\[controller\]=DemoLicensesGeneration&cHash=3f31f0da1fc6e9f7244fc60490596194](https://www.bruker.com/service/support-upgrades/software-downloads/nmr/free-topspin-processing-...olicensesgeneration[controller]=DemoLicensesGeneration&cHash=3f31f0da1fc6e9f7244fc60490596194)

## Step 4: Now, you should write down or copy your “Personal license code” for later usage.

Free NMR Topspin License P

https://www.bruker.com/service/support-upgrades/software-downloads

BRUKER

Products Applications Service News Events About Us Logout Contact

Free TopSpin Processing for Academia

Contact

Home . Service . Support & Upgrades . Software Downloads . NMR . NMR TopSpin License for Academia

### NMR TopSpin License Processing for Academia

In order to provide students, researchers and teachers with unlimited access to the best tools for off-line NMR processing, Bruker is making their market leading NMR processing software TopSpin available free of charge for all academic users. In addition to Bruker NMR data, the software is able to process data acquired on Agilent, Varian and Jeol systems, as well as read any data provided in JCAMP format. We are happy to offer you a free Academia license for TopSpin 4 and TopSpin 3.6 ("Academia license").

#### Free License For Academia

With the following code you can activate your free academia license:

**4WBSB-XXXX-XR8DT-XXXX-D4VKZ**

1. Write down this code. (Personal license code)

Just write it down (or copy it to the clipboard), start TopSpin, choose the option "Activate this installation now" and follow the instructions. If you have not installed TopSpin already you can download it [here](#).

Your license is valid for 15 years. You can use it on a compatible computer of your choice and you can move it to another system if needed. See the section "Rehosting of license" in the license management manual for details.

[Download for Windows](#)   [Download for Linux](#)   [Download for Mac OS](#)

2. Click here to download software for Windows.

Disclaimer

Access to Academia licenses is restricted to academic customers. By using this web page for license requests you confirm that you are employed by or

# Step 5: Download & Install TopSpin & CMC-assist. (select “Data processing only” when installation)

**TopSpin - Windows PC**

**Current versions:**

- TopSpin 4.0.7 (for AVANCE NEO and processing in general)
- TopSpin 3.6.1 (for AVANCE III, III HD and processing in general)
- TopSpin 3.5 pl7 (for AVANCE III, III HD and processing in general)
- TopSpin 3.2 pl7 (for AVANCE II, III, III HD, Fourier 300 and processing in general)
- TopSpin 2.1 pl8 (for AVANCE I)
- TopSpin 1.3 pl10 (for AVANCE DRX, DPX, DMX, and DSX)

For detailed information refer to the corresponding [release letter](#) and [software support matrix](#).

(For information on acquiring the current version of TopSpin on DVD, contact your nearest [Bruker BioSpin office](#).)

To work with TopSpin a license is required. To download the current TopSpin version, click on the link and follow the instructions provided below. Please store the file on your hard disk and start the installation when the download is complete.

**Current Release Download**

**TopSpin 4.0.7 & CMC-assist 2.14 (1718.9 MiB)**  
(Created: 2019-05-07)  
TopSpin for AVANCE NEO and processing in general.  
This link allows for the direct installation of a complete acquisition/processing version for Windows 7 and Windows 10.  
MD5 checksum:  
ed3bb65b45ac8759970e481a6a46d75d

**TopSpin 3.6.1 & CMC-assist 2.13 (990.2 MiB)**  
(Created: 2018-12-21)  
TopSpin for AVANCE III, III HD and processing in general.  
This link allows for the direct installation of a complete acquisition/processing version for Windows 7 and Windows 10.  
Very large self-extractable file --> fast connection required!

**TopSpin 3.2 pl7 & CMC-assist 2.1 (732.1 MiB)**  
(Created: 2015-07-21)  
TopSpin for AVANCE II, III, III HD, Fourier 300 and processing in general.  
This link allows for the direct installation of a complete acquisition/processing version for Windows 7

**ISO - Downloads**

These images contain everything that you would also see on a DVD when you order it from Bruker. They contain all programs for all supported operating systems and also all additional software like AMIX, AURELIA and CMC-assist. For programs that are supported on the MAC, the MAC version is also contained in this image.

**TopSpin 4.0.7 (6441.5 MiB)**  
(Created: 2019-05-07)  
For Avance NEO.  
MD5 checksum:  
f74f4220a82fc427eae5c1f674684e3


**TopSpin 3.6.1 (rel36.iso) (4259.3 MiB)**  
(Created: 2018-12-21)  
For Avance III, III HD.  
MD5 checksum:  
6811dac59a729ddf4d5bda97dd609b11

**TopSpin 3.2 pl.7 (rel32.iso) (3855.4 MiB)**  
(Created: 2015-07-21)  
For Avance II, III, III HD, Fourier 300

[↑](#)  
[f](#)  
[t](#)  
[in](#)  
[✉](#)

[https://www.bruker.com/nc/service/support-upgrades/software-downloads/nmr/pc-topspin.html?...ercontent\\_jwtfiledownload\[controller\]=JwtFileDownload&cHash=d2e1cd7b2c97b53c864ef9ba21ea6b4](https://www.bruker.com/nc/service/support-upgrades/software-downloads/nmr/pc-topspin.html?...ercontent_jwtfiledownload[controller]=JwtFileDownload&cHash=d2e1cd7b2c97b53c864ef9ba21ea6b4)

## Step 6: After TopSpin installation, start TopSpin and install License ticket.

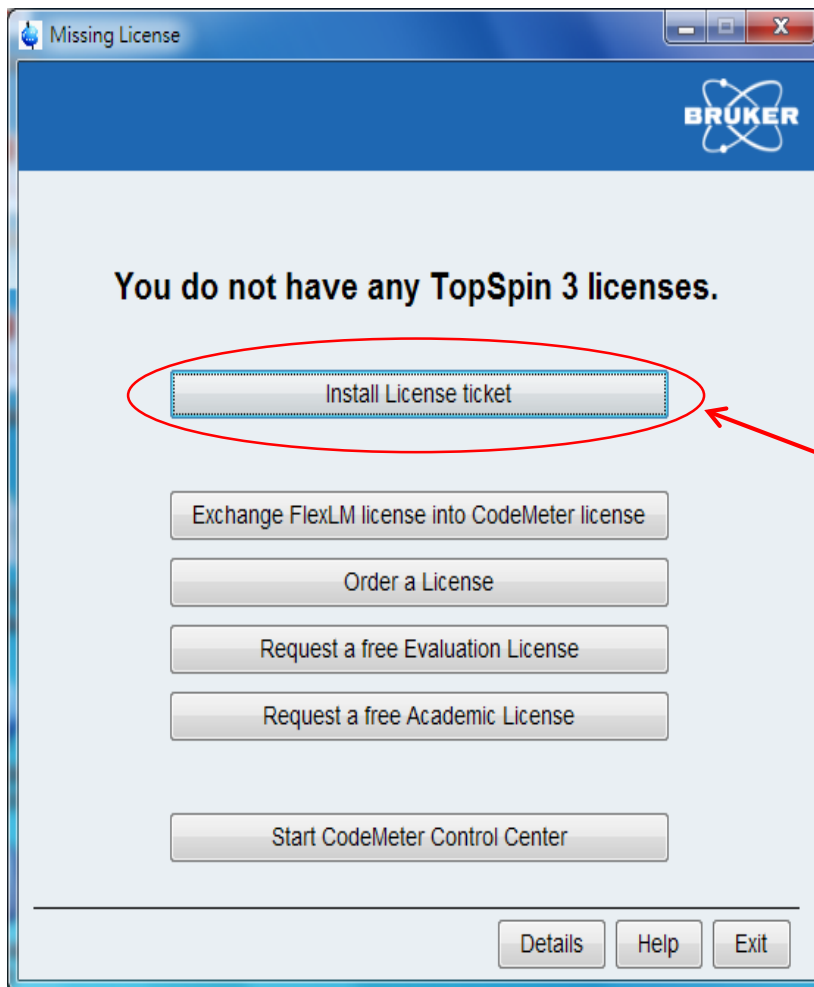


Click "Install License ticket".


NOTE: Virus protection software may cause trouble during installation, you may need to stop the protection for a while.

## Step 6.1

You should see the welcome screen of the activation process in browser :

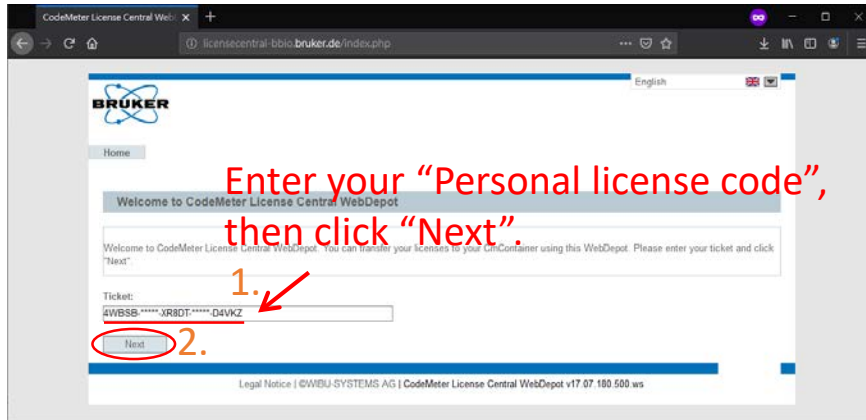


## Step 6.2

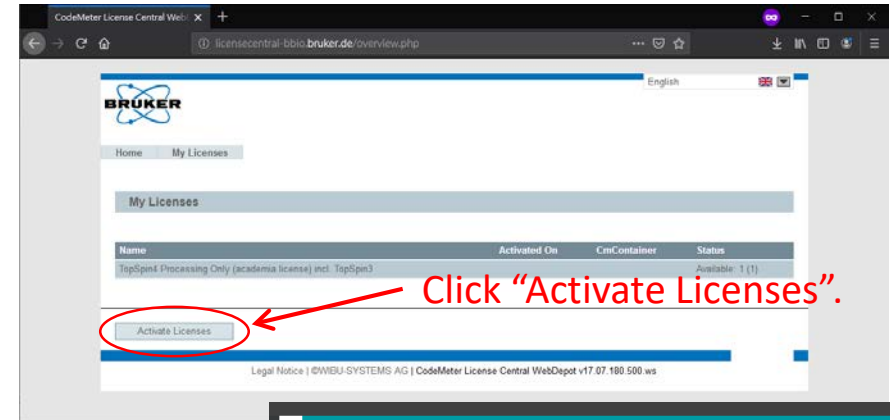

## Step 6.3

You should select the right option "CmActLicense"

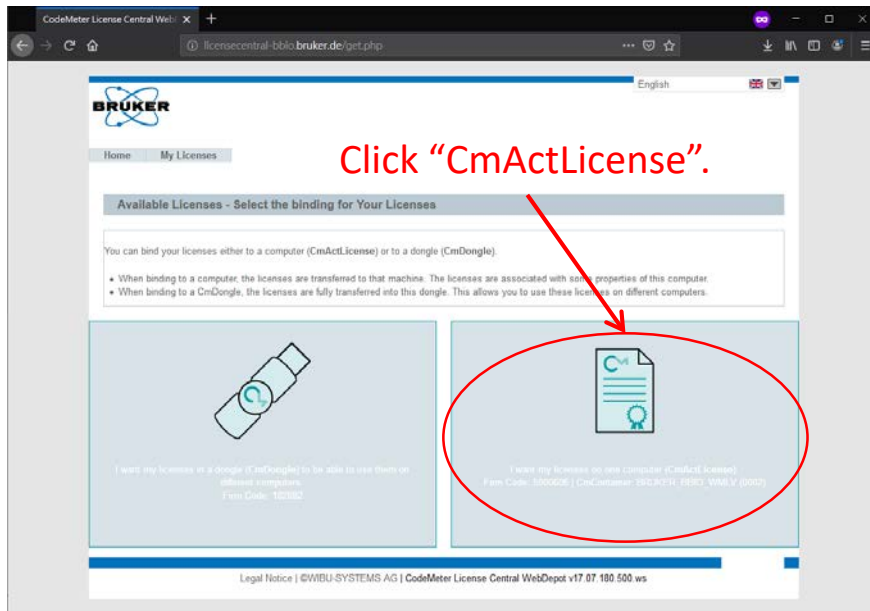

## Step 6.4

