# Free Topspin & CMC Installation Guide 學術版免費 Topspin & CMC 安裝步驟 2018.05.25

If you already have free Topspin3.5 pl7 with CMC\_assist2.6, you don't need to reinstall it (but will expire on 2019.03.31).

Step 1: Register for Bruker users account (需先註冊才能取的免費license)

https://www.bruker.com/about-us/register.html

RUKER	Products Applications	Service News Events About us	Lopin Q	
Home - Aboutus - R	itur.			
Register Welcome to Bruke software downloa	r for Bruker user r quick registration portal. To access t d, you are required to have an accoun	S account o some of our Bruker community information such a	s user documentation, events, newsletter,	
if you don't have a	n account with us yet, you can start n	ow by filling out the form below.		
First Name: *		I am interested in the following n  Magnetic Resonance eNews i  Midanular Vibrational Source	evaletters: . The Resonance Newscast	
Country or Area: *	Talwan	Mass Spectrometry		
Email Address: * Create Password: *		FIRST Newsletter (X-ay Diffra     Nanovations (Atomic Force N	ction and Elemental Analysis) (croscopy)	
Confirm Password		Surface Insights (Surface Pro	(ling and Mechanical Testing)	
Password must be	at least 6 characters long	Preclinical Imaging energy		
		BioViews (Life Sciences Micro 	scopy)	
		CBRNE Detection e-nexis		
		Microbiology & Diagnostics		

### Step 2: Login using your account , and click "Free Academia license" (至下面網頁, 登入後 點選 Free Academia license)

https://www.bruker.com/service/support-upgrades/software-downloads/nmr.html

Home + Senice + Support & Upgrades + Software Downloa	ás + NNR	
NMR Software & Dov	vnloads	🖌 Log in
This section is available for customer only, registe	red users can access to NMR software download, upgra	dy, and patches for Windows, Linux and MAC OS X.
If you don't have an account, please register for a listed software products below.	n account, if you already have an account pease log in h	are to access the download section containing all
Software Downloads	Generate License Code	License Request
If you are logged in, you can download the software below by clicking on the link:	TopSpin processing software is free for academia and government institutions. And you	Request NMR processing and acquisition licenses
<ul> <li>TopSpin for Windows, Linux and MAC OSX TopSpin installation includes following products:</li> </ul>	can get an Evaluation license for the full Bruker. Suite here (requires login/ quick registration).	Request Form
iconNMR	Free Evaluation license	Choose "Free
SmartDriveNMR     CMC a		Academia license"
CMC-se		
CMC assist		
Potencysee     FBS		
TopSolids		
<ul> <li>BioTop</li> </ul>		
<ul> <li>APSY</li> </ul>		
<ul> <li>NMR-SIM</li> </ul>		
<ul> <li>AssureNMR for Windows and Linux</li> </ul>		

## Step 3: Now, Request "Personal license Ticket" (進入後, 點選" Request personal license ticket now")



# Step 4: Now, you should write down or copy your "Personal license code" for later usage (記下你的專屬" license code", 之後會用到)



### Step 5: Download & Install (process only) Topsoin & CMC 下載 Topspin &, 並照步驟安裝 (process 功能即可)

Free TopSpin Processing for Acader	nia	TopSpin - Windows PC	
	Provide State	Current versions:	(For information on acquiring the current version of TopSpin on DVD, control of a second Top Spin Price of Sci
Contact rares - Berlin - Bartert Scapada - Bathwert Scientiada - Villi - Par Tatlor	nowny - 107 Tatler Canada Frances	TopDari 4.0.2 (WHAVECE NED and processing in general     TopDari 4.0.2 (WHAVECE NED and processing in general     TopDari 1.2 or // WHAVECE (W, WHAVECE W, W, WHAVECE W, W, WHAVECE W,	sumation memory on the second se
NMR Topspin License Process	ing for Academia	Current Release Download	
In order to provide students, researchers and teachers with unlimited a leading MMR processing software TooSpin available free of charge for a data acquired on Aglient, Varien and Jeol systems, as well as read any	coss to the best tools for off-line NMR processing, Bruker is making their marker, I academic users. In addition to Bruker MMR data, the otherwe is able to process ats provided in JCAMP format, We are happy to offer you a free Academia Ioanse for	Toptor 4036 CVC-west 210 (493.7 MB)	ISO - Downloads
TopSpin 4 ('Academia license').		(Created: 2018-05-15) TooSpin for XIANCE NED and processing in general	These images contain everything that you would also see on a DiD when you order it from Bruker. They contain all programs for all succorted
Free License For Academia		This link allows for the direct installation of a complete acquisition/processing version for Windows 7 and Windows 10.	operating systems and also all additional software like AMU, AUREUA and CMC-assist. For programs that are supported on the MAC, the MAC version
With the following code you can activate your free academia license		MDS checksum P1x95666097cb1483566646555ba3ec2	is also contained in this image.
HNL56	BUV4S-EN3Y6		TopSpin 4.0.3 (mi40.iso) (6146.9 Mi8)
Just write it down (or copy it to the cipboard), start ropopin, choose to installed Top5pin already you can download it here.	eloption individual this insclusioon now "and toxice the instructions, if you have not	Detain 350100Cause 25(966.6 MB)	(Created: 2018-03-15) For Assess NEC
Your license is valid for 15 years. You can use it on a compatible compu- "Rehoting of license" in the license management manual for details.	ar of your choice and you can move it to another system if needed. See the section 🔹 🚹	(Created 2016-05-12) TopDon for ANANCE II., II HD and processing in general. This link allows for the direct installation of a complete	NDS checkaum: 4117dx1e54666644245a95e48429926
Download for Windows Download for Lin	un Downlaad for Nac OS	acque don processing version to mindows 2 and mindows 20. Very large self-extractable file —> fast connection required	TopSpin 3.6.0 (w/34.0d) 4410.2 Mills
Fyou need the Academia version of TopSpin 3.5 instead, please follow	the line in	27a38cH4ca7852810c1424815a96c199	(Created: 2018-03-13)

Topspin3.6 with free NUS processing . New license code will valid for 15 years. Step6: After install topspin package, Click your topspin icon to install license ticket 安裝後, 第一次開啟, 會出現以下畫面, 點選 install license ticket, 然後照步驟完成 license

Missing Licens	e BRUKEF	
You	do not have any TopSpin 3 licenses.	Click to activate your free license
	Exchange FlexLM license into CodeMeter license	
	Request a free Evaluation License Request a free Academic License	
	Start CodeMeter Control Center	
	Details Help Exit	

NOTE: Virus protection software may cause trouble during installation, you may need to stop the protection for a while 某些防毒軟體, 或網頁安全設定可能會阻止安裝, 請先暫停防毒軟體及允許網頁安全

#### Step6.1

Ston6 3	
Stepo.5	Select the binding for Your Licenses



### Step6.2

The activation services will now display the license package(s) which have been deposited for the provided ticket.

Address   passa   assis (Bellance al 1.18 107 Mit - Electron Internet   ad one		- (D - A
🕞 🖉 trg//sensental this balance includes the set dependence of the	• 4 X 3	ing P
norther 😰 CodeMater Lineaus Central WebDeps		
$\sim$		
UKER		
me Myticaniel	English	90 E
My Licenses		
uni	Solution	Talat
RUNZER FLERONLEV Single Loursequede Andreits (and in Wideys)		Anilde
Attentioner (6.2) Click "Act	ivata Liconso"	
(0.2) CIICK AC		
Legal Rotor ( & WEG SYSTEMS AD / CoduMeter Loanse Central We	BCxpct v15.08.152.508 (2016-01-20.10.38.48 (J/TC)	

You should select the right option "CmActLicense".

